
InPro 6850 i, 6900 i & 6950 i Series O₂ Sensors

Instruction manual
Bedienungsanleitung
Instructions d'utilisation

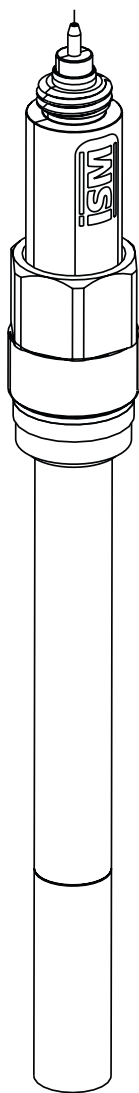InPro 6850i/6900i/6950i
52 206 349 G

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InPro 6850 i, 6900 i & 6950 i Series O₂ Sensors

Instruction manual

InPro, InDip, InFit, ISM and InTrac are trademarks
of the METTLER TOLEDO Group.

Contents

| | | |
|------------|---|-----------|
| 1. | Introduction | 5 |
| 2. | Important notes | 6 |
| 2.1 | Notes on operating instructions | 6 |
| 2.2 | Intended use | 6 |
| 2.3 | Safety instructions | 7 |
| 2.4 | Environmental protection | 8 |
| 2.5 | Examples of some typical applications | 9 |
| 2.6 | Use in Ex-zones | 9 |
| 2.7 | Ex-classification ATEX | 10 |
| 2.7.1 | Introduction | 10 |
| 2.7.2 | Rated data | 10 |
| 2.7.3 | Special conditions | 11 |
| 2.8 | Ex-classification FM approved | 12 |
| 2.8.1 | Introduction | 12 |
| 2.8.2 | Rated data | 13 |
| 2.8.3 | Special conditions | 13 |
| 2.8.4 | Applied standards: | 14 |
| 3. | Product description | 15 |
| 3.1 | General information | 15 |
| 3.2 | Principle | 15 |
| 3.3 | Scope of delivery | 16 |
| 3.4 | Equipment features | 17 |
| 4. | Installation | 18 |
| 4.1 | Mounting the sensor | 18 |
| 4.2 | Connection | 18 |
| 4.2.1 | Connecting the sensor to an AK9 cable | 18 |
| 4.2.2 | Connecting the AK9 cable to the transmitter | 19 |
| 5. | Operation | 20 |
| 5.1 | Start-up and polarizing | 20 |
| 5.2 | Calibration | 21 |
| 5.2.1 | Purpose of calibration | 21 |
| 5.2.2 | What you have to know for calibration | 21 |
| 5.2.3 | Single point calibration | 22 |
| 5.2.4 | Dual point calibration | 22 |
| 6. | Maintenance | 23 |
| 6.1 | Inspection of the sensor | 23 |
| 6.1.1 | Visual inspection | 23 |
| 6.1.2 | Testing the METTLER TOLEDO O ₂ sensor master | 24 |
| 6.1.3 | Testing the sensor via a transmitter | 25 |
| 6.1.4 | ISM design | 26 |
| 6.2 | Changing the electrolyte, the membrane body or the interior body | 26 |
| 7. | Storage | 29 |
| 8. | Product specification | 29 |
| 8.1 | Certificates | 29 |
| 8.2 | Specifications | 30 |
| 9. | Ordering information | 32 |
| 9.1 | Sensors with ISM functionality | 32 |
| 9.2 | Accessories | 32 |
| 9.3 | Spare parts | 32 |
| 9.4 | Recommended transmitters | 33 |
| 9.5 | Recommended housings | 34 |
| 10. | Theory of the polarographic sensor | 35 |
| 10.1 | Introduction | 35 |
| 10.2 | Principle of the design of an oxygen electrode | 35 |
| 10.3 | Parameters determining current | 39 |
| 10.4 | Polarization voltage | 39 |
| 10.5 | Temperature | 40 |
| 10.6 | Dependence on flow | 40 |
| 10.7 | Oxygen partial pressure – oxygen concentration | 40 |

1 Introduction

Thank you for buying this **InPro™ 6850 i/6900 i/6950 i sensor from METTLER TOLEDO.**

The construction of the InPro series employs leading edge technology and complies with safety regulations currently in force. Notwithstanding this, improper use could lead to hazards for the user or a third-party, and/or adverse effects on the plant or other equipment.

Therefore, the operating instructions must be read and understood by the persons involved before work is started with the sensor.

This instruction manual must always be stored close at hand, in a place accessible to all people working with the InPro sensor.

If you have questions, which are not or insufficiently answered in this instruction manual, please contact your METTLER TOLEDO supplier. He will be glad to assist you.

2 Important notes

2.1 Notes on operating instructions

These operating instructions contain all the information needed for safe and proper use of the InPro 6850i / 6900i / 6950i sensor.

The operating instructions are intended for personnel entrusted with the operation and maintenance of the sensor. It is assumed that these persons are familiar with the equipment in which the sensor is installed.

Warning notices and symbols

This instruction manual identifies safety instructions and additional information by means of the following symbols:

This symbol draws attention to safety instructions and warnings of potential danger which, if neglected, could result in injury to persons and/or damage to property.

This symbol identifies additional information and instructions which, if neglected, could lead to defects, inefficient operation and possible loss of production.

2.2 Intended use

METTLER TOLEDO InPro 6850i/6900i/6950i sensors are intended solely for inline measurement of the oxygen partial pressure in liquids and gases, as described in this instruction manual.

Any use of these sensors which differs from or exceeds the scope of use described in this instruction manual will be regarded as inappropriate and incompatible with the intended purpose.

The manufacturer/supplier accepts no responsibility whatsoever for any damage resulting from such improper use. The risk is borne entirely by the user/operator.

Other prerequisites for appropriate use include:

- compliance with the instructions, notes and requirements set out in this instruction manual.
- acceptance of responsibility for regular inspection, maintenance and functional testing of all associated components, also including compliance with local operational and plant safety regulations.
- compliance with all information and warnings given in the documentation relating to the products used in conjunction with the sensor (housings, transmitters, etc).
- observance of all safety regulations governing the equipment in which the sensor is installed.
- correct equipment operation in conformance with

the prescribed environmental and operational conditions, and admissible installation positions.

- consultation with METTLER TOLEDO Process Analytics in the event of any uncertainties.

2.3 Safety instructions

- The plant operator must be fully aware of the potential risks and hazards attached to operation of the particular process or plant. The operator is responsible for correct training of the workforce, for signs and markings indicating sources of possible danger, and for the selection of appropriate, state-of-the-art instrumentation.
- It is essential that personnel involved in the commissioning, operation or maintenance of these sensors or of any of the associated equipment (e.g. housings, transmitters, etc.) be properly trained in the process itself, as well as in the use and handling of the associated equipment. This includes having read and understood this instruction manual.
- The safety of personnel as well as of the plant itself is ultimately the responsibility of the plant operator. This applies in particular in the case of plants operating in hazardous zones.
- The oxygen sensors and associated components have no effect on the process itself and cannot influence it in the sense of any form of control system.
- Maintenance and service intervals and schedules depend on the application conditions, composition of the sample media, plant equipment and significance of the safety control features of the measuring system. Processes vary considerably, so that schedules, where such are specified, can only be regarded as tentative and must in any case be individually established and verified by the plant operator.
- Where specific safeguards such as locks, labels, or redundant measuring systems are necessary, these must be provided by the plant operator.
- A defective sensor must neither be installed nor put into service.
- Only maintenance work described in this operating instruction may be performed on the sensors.
- When changing faulty components, use only original spare parts obtainable from your METTLER TOLEDO supplier (see spare parts list, "Section 9.3").
- No modifications to the sensors and the accessories are allowed. The manufacturer accepts no responsibility for damages caused by unauthorized modifications. The risk is borne entirely by the user.

- When using cables for sensor connection not supplied and recommended by the manufacturer the following standards need to be considered:
 - US National Electrical Code® (ANSI/NFPA 70 [NEC®]), where applicable.
 - Canadian Electrical (CE) Code® (CEC Part 1, CAN/CSA-C22.1), where applicable.
- For guidance on US installations, see ANSI/ISA-RP12.06.01, Installation of Intrinsically Safe Systems for Hazardous (Classified) Locations.
- Control room equipment connected to intrinsically safe associated apparatus in the US should not use or generate more than the specified Um of the associated apparatus.
- Care must be taken during installation to avoid impacts or friction that could create an ignition source.
- Tampering and replacement with non-factory components may adversely affect the safe use of the system.
- Insertion or withdrawal of removable electrical connectors or modules is to be accomplished only when the area is known to be free of flammable vapors.
- **WARNING** –
INTRINSICALLY SAFE APPARATUS CAN BE A SOURCE OF IGNITION IF INTERNAL SPACINGS ARE SHORTED OR CONNECTIONS OPENED.
- **WARNING** –
DO NOT OPEN WHEN AN EXPLOSIVE ATMOSPHERE IS PRESENT.
- **WARNING** –
SUBSTITUTION OF COMPONENTS MAY IMPAIR INTRINSIC SAFETY.
- **WARNING** –
SUBSTITUTION OF COMPONENTS MAY IMPAIR SUITABILITY OF THE EQUIPMENT
- **WARNING** –
FOR CONNECTION ONLY TO NON-FLAMMABLE PROCESSES.
- **WARNING** –
TO MAINTAIN THE ENCLOSURE IP66 RATING, THE CONNECTOR MUST BE FULLY ENGAGED.

2.4 Environmental protection

Waste electrical products should not be disposed of with household waste. Please recycle where facilities exist. Check with your Local Authority or retailer for recycling advice.

2.5 Examples of some typical applications

Below is a list of examples of typical fields of application for the oxygen sensors. This list is not exhaustive.

Measurement in liquids:

- Biotech
- Chemical applications
- Brewing
- Beverage filtration
- Filling stations

Measurement in gases:

- CO₂ recovery
- CO₂ purity
- Product storage
- Inert production

2.6 Use in Ex-zones

Attention! For an installation in Ex-zones please read the guidelines following hereafter:

Ex-classification ATEX:

⊠ II 1/2G Ex ia IIC T6...T3 Ga/Gb

⊠ II 1/2D Ex ia IIIC T69 °C...T161 °C Da/Db

Number of the test certificate:

SEV 14 ATEX 0169X

IECEX SEV 14.0026X

CML 22 UKEX 2212X

Ex-classification FM approved:

IS/I, II, III/1/ABCDEFG/T6 Ta = 60 °C
– 53 800 002; Entity

Project ID: 3021227

FM certificate number: FM16US0034X

FM18CA0021X

2.7 Ex-classification ATEX

2.7.1 Introduction

- According to Directive 2014/34/EU (ATEX 114)¹⁾ Appendix I, InPro 6XXX*/**/* oxygen sensors are devices group II, category 1/2G and according to RL 99/92/EG (ATEX 137)²⁾ may be used in zones 0/1 or 1/2 or 1 or 2 and gas groups IIA, IIB and IIC that are potentially explosive due to combustible substances in above stated temperature classes.
- For use / installation, the requirements of EN 60079-14 must be observed.
- According to Directive 2014/34/EU (ATEX 114)¹⁾ Appendix I, InPro 6XXX*/**/* oxygen sensors are devices group II, category 1/2D and according to RL 99/92/EG (ATEX 137)²⁾ may also be used in zones 20/21 or 21 or 22 that contain combustible dusts.
- The digital sensor circuit is part of a common intrinsically safe system and is for operation connected to a separately certified transmitter.
- The digital sensor circuit as part of an intrinsically safe system is isolated from the not-intrinsically safe electric circuits up to a maximum rated voltage of 375 V and from grounded parts up to a maximum rated voltage of 30 V.

2.7.2 Rated data

Measuring circuit

Intrinsically safe ignition protection class Ex ia IIC only for connection to a certified intrinsically safe circuit.

Digital O₂ oxygen sensor maximum values

Two-wire current circuit:

$$U_i \leq 16 \text{ V}$$

$$I_i \leq 30 \text{ mA}$$

$$P_i \leq 50 \text{ mW}$$

$$L_i = \text{negligible}$$

$$C_i = \text{negligible}$$

¹⁾ For UK statutory Requirements SI 2016 No. 1107

²⁾ For UK regulations 7 and 11 of the Dangerous Substances and Explosive Atmospheres Regulations 2002 (DSEAR)

2.7.3 Special conditions

- The relationship between the maximum permissible ambient or media temperature and temperature class, for category 1G applications, zone 0, is shown in the following table:

| Temperature class | Max. ambient or media temperature |
|--------------------------|--|
| T 6 | 68 °C |
| T 5 | 80 °C |
| T 4 | 108 °C |
| T 3 | 160 °C |

- The relationship between the maximum permissible ambient or media temperature and temperature class, for category 1D applications, zone 20, is shown in the following table:

| Temperature class | Max. ambient or media temperature |
|--------------------------|--|
| T 69 °C | 68 °C |
| T 81 °C | 80 °C |
| T 109 °C | 108 °C |
| T 161 °C | 160 °C |

- The capacitance and inductance of the connecting cable has to be considered.
- The O₂ oxygen sensor type InPro 6XXX can be used in/with the fittings InFit™ 76*_*_*_* or InTrac™ 7**_*_*_*, or in/with other suitable fittings in potentially explosive areas.
- The metal body of the O₂ oxygen sensors, or the fittings InFit 76*_*_*_* or InTrac 7**_*_*_*, or other appropriate fitting is optionally included in the routine pressure test of the system.
- The independent fitting used for installation of the O₂ oxygen sensor must be conductively connected to the equipotential bonding system.


2.8

Ex-classification FM approved

2.8.1 Introduction

The following FM control drawing and the standards listed in section 2.8.4 must be observed, where applicable: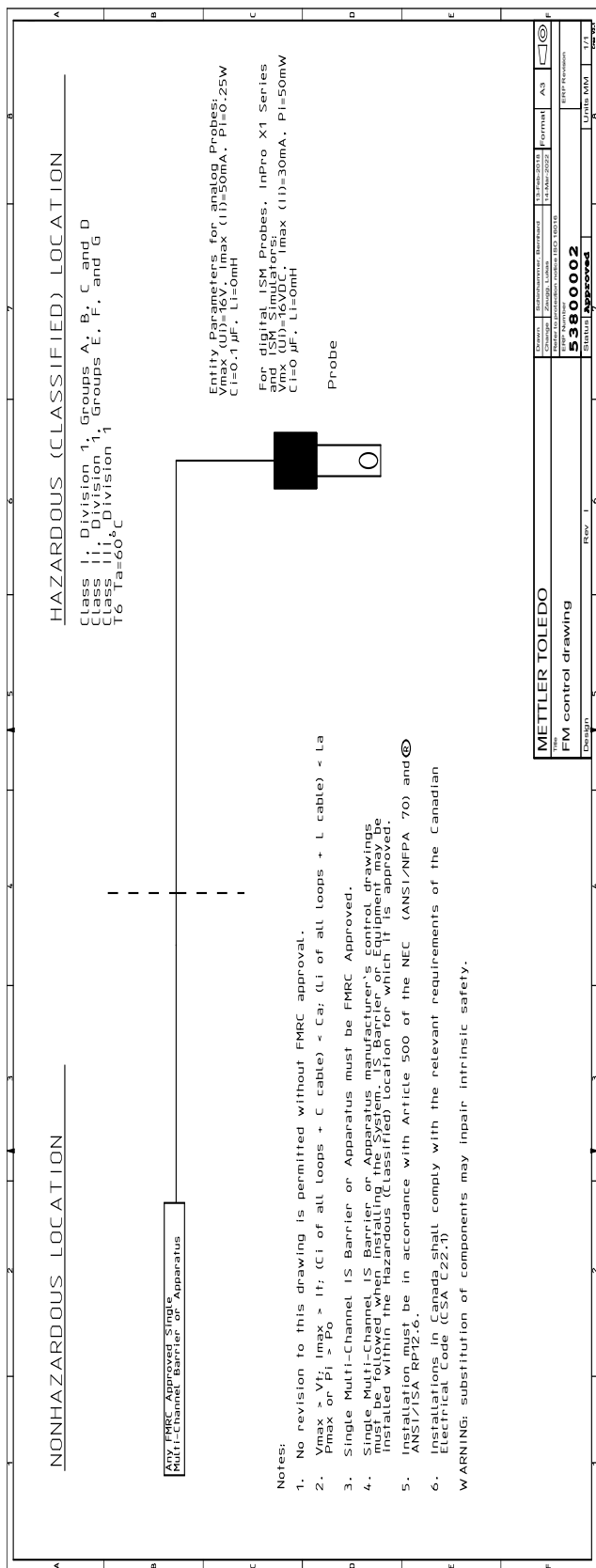

2.8.2 Rated data

For rated data please observe section 2.7.2

2.8.3 Special conditions

In intrinsically safe apparatus, the analytical probes (InPro 6000 Series with InFit 76X/Y and InTrac 7XX/YY Series housings) equipment is designated with the following specific conditions of use.

1. The oxygen sensor shall be installed in compliance with the enclosure, mounting, spacing and segregation requirements of the ultimate application, including a tool removable cover, and is suitable for use with FM Approved InFit 76X/Y and InTrac 7XX/YY Series housings.
2. Process temperature no greater than +130 °C.
3. Maximum permissible working pressure is 12 barg (174 psig).
4. Potential Electrostatic Charging Hazard – To prevent the risk of electrostatic sparking, the non-metallic surface should only be cleaned with a damp cloth.
5. Enclosures containing titanium constitute a potential risk of ignition by impact or friction. Care must be taken during installation and use to prevent impact or friction.

2.8.4 Applied standards:

United States Standards

| Title | Number | Issue Date |
|--|-----------------------|-------------------|
| Approval Standard for Electrical Equipment for Use in Hazardous (Classified) Locations – General Requirements | FM Class 3600 | 2022 |
| Approval Standard for Intrinsically Safe Apparatus and Associated Apparatus for Use in Class I, II & III, Division 1, Hazardous (Classified) Locations | FM Class 3610 | 2021 |
| Approval Standard for Electrical Equipment for Measurement, Control and Laboratory Use | FM Class 3810 | 2021 |
| Explosive Atmospheres – Part 0: Equipment – General Requirements | ANSI/UL 60079-0:2020 | 2020 |
| Explosive Atmospheres – Part 11: Equipment Protection by Intrinsic Safety “i” | ANSI/UL 60079-11:2018 | 2018 |
| Safety Requirements for Electrical Equipment for Measurement, Control, and Laboratory Use – Part 1: General Requirements | ANSI/UL 61010-1:2018 | 2018 |

Canadian Standards

| Title | Number | Issue Date |
|--|-------------------------------------|-------------------|
| Explosive Atmospheres – Part 0: Equipment – General Requirements | CAN/CSA-C22.2 No. 60079-0:2019 | 2019 |
| Explosive Atmospheres – Part 11: Equipment Protection by Intrinsic Safety “i” | CAN/CSA-C22.2 No. 60079-11-14:R2018 | R2018 |
| Safety Requirements for Electrical Equipment for Measurement, Control, and Laboratory Use – Part 1: General Requirements | CAN/CSA-C22.2 No. 61010-1-12:R2017 | R2017 |

3 Product description

3.1 General information

The **oxygen sensor series InPro 6850 i/6900 i/6950 i** with integrated temperature probe are used for measurement of oxygen at low and medium concentrations.

The sensors are **sterilizable**, and in the majority of cases are also autoclavable. They are **compatible with CIP** (Cleaning-In-Place).

InPro 6xxx i sensors with ISM™ functionality offer Plug and Measure as well as enhanced diagnostics features.

3.2 Principle

Amperometric oxygen sensors:

The amperometric oxygen sensors of METTLER TOLEDO are all based on the Clark measuring principle. However different series are offered, which differ in the number and arrangement of their electrodes and thus in their specifications:

- The measurement system of the InPro sensors consists of a working electrode (cathode), a counter electrode (anode), a reference electrode and a guard ring electrode. The measurement system is separated from the process medium by an oxygen permeable membrane.
- The measuring technique in principle is alike with all sensors.
- The transmitter supplies a constant voltage between cathode and anode.
- The electrolyte creates a conductive connection between the electrodes.
- The oxygen molecules migrate from the measurement medium through the membrane to the cathode to which the voltage is supplied and are reduced. At the same time, oxidation takes place at the anode.
- This causes a current to flow between the anode and cathode which is directly proportional to the partial pressure of oxygen (pO₂) in the process medium. With digital ISM™ sensors, the sensor converts the current into oxygen concentration and communicates the value to the transmitter.
- The guard ring electrode of the InPro 6900 i and 6950 i built into the sensor reduces the oxygen that migrates to the side of the cathode and that can distort the measurement. The guard ring electrode therefore enables traces of oxygen to be precisely determined at even the lowest concentrations.

Note: Please refer to “Section 10 – Theory of the polarographic sensor” for further information.

ISM sensors:

All oxygen sensors with the index "i" (6850i, 6900i, 6950i) are equipped with ISM.

Principle: In the sensor head a chip is integrated, which takes over the entire monitoring and control of the sensor and, beyond that, stores all sensor data.

The following data are available permanently in the sensor:

- type of sensor
- serial no.
- software version
- hardware version
- order no.
- operating time
- calibration time and calibration date
- calibration table

To check the system, the following indicators are supervised:

- temperature
- slope
- zero current
- air current
- polarization voltage

Based on this information, the transmitter calculates a wear monitor and displays it depending upon transmitter type. (see the respective instruction manuals)

ISM enables the connection of the digital sensor to iSense. This user-friendly software allows management of all sensor information and storage of it in a database. Furthermore, the sensor can be calibrated via the iSense software.

3.3 Scope of delivery


Each sensor is supplied fully assembled and factory-tested for correct function together with:

- a quality control certificate
- inspection certificates 3.1
(complying with EN 10204)

METTLER TOLEDO ISM O₂ sensors are supplied with fitted membrane body but without electrolyte, covered with the protection cap. They have been checked for proper functionality. See electrolyte order details at section "Spare Parts", 9.3.

3.4 Equipment features

12 mm sensor

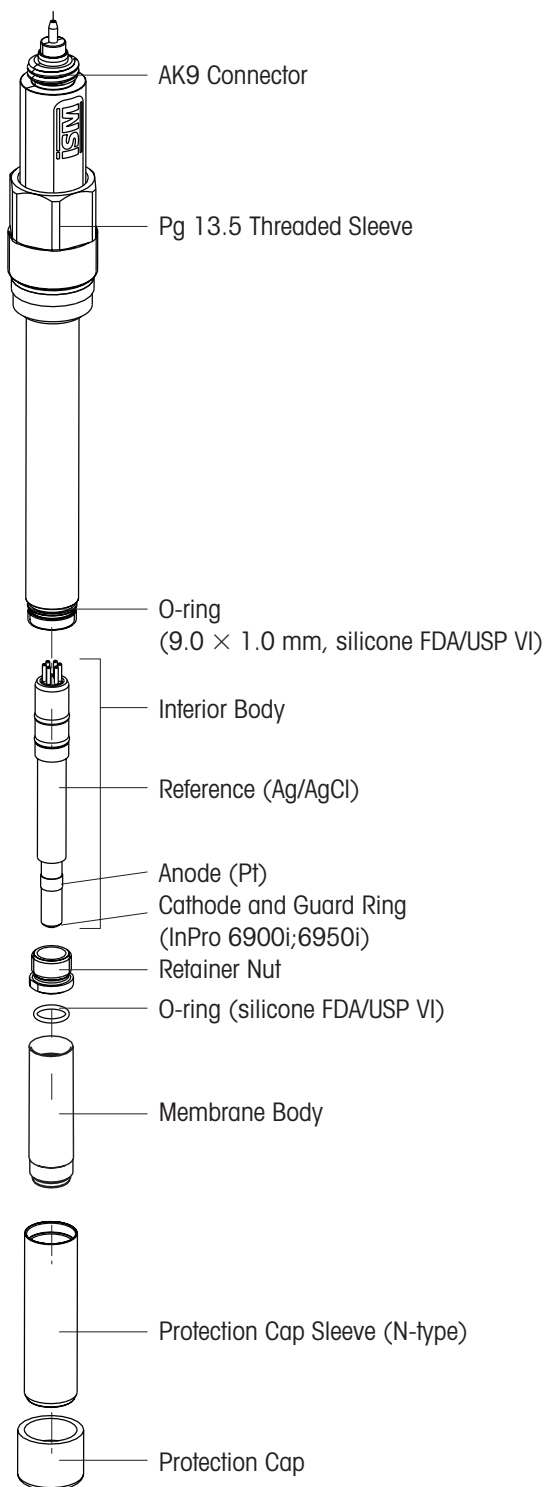4 Installation

4.1 Mounting the sensor


Important! Remove the protection cap before mounting the sensor.

Mounting the sensor in a housing

Please refer to the instruction manual of your housing explaining on how to mount the sensor in place.

Mounting the sensor directly in a pipe or a vessel

The 12 mm sensors can be mounted directly through a socket with inside thread Pg 13.5 and securely tightened via the Pg 13.5 threaded sleeve.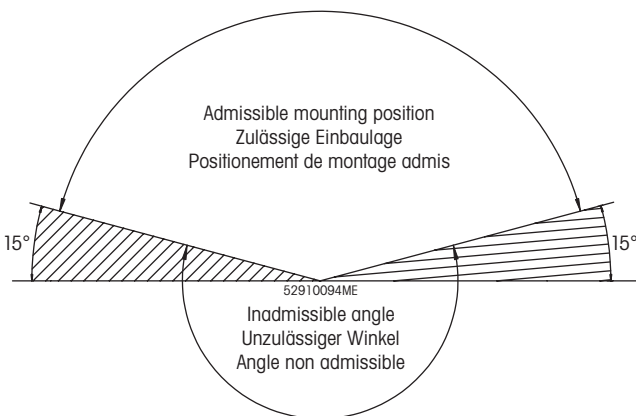

4.2 Connection

4.2.1 Connecting the sensor to an AK9 cable


The sensor is connected to the transmitter via an AK9 cable. The AK9 cable ensures a secure connection between the transmitter and the sensor under harsh industrial conditions. The robust watertight IP 68 connector housing ensures maximum process safety.


Do not touch the sensor at the AK9 connector plug!

Tightly screw the plug to fasten the two parts.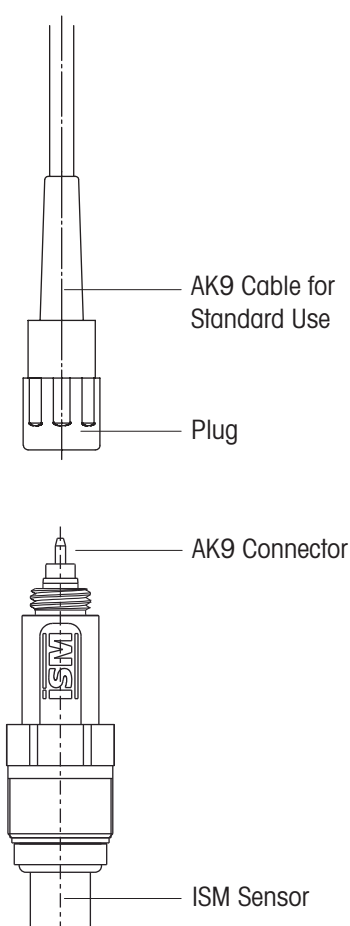

4.2.2 Connecting the AK9 cable to the transmitter

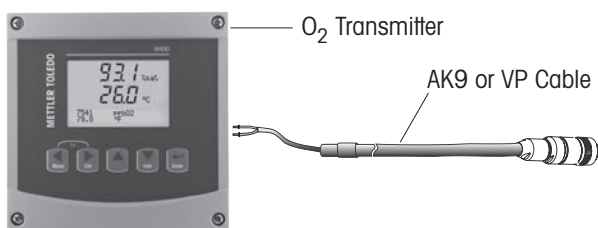Note: For connecting the AK9 cable to the terminals of the transmitter, please refer to the instructions given in the METTLER TOLEDO transmitter manual.

5 Operation

Important! Before using the sensor for the first time, electrolyte must be added to the membrane body (see “Chapter 6.2”).

5.1 Start-up and polarizing

Important! The protection cap must be removed before mounting the sensor in the process.

When the system is operated for the first time or if the sensor has been disconnected from the voltage source (transmitter or O₂ sensor master) for longer than 5 minutes, the sensor has to be polarized prior to calibration by connecting it to the operating O₂ transmitter or to a sensor master. After 6 hours, the sensor is fully polarized and ready for operation.

Attention: The InPro 6950 i should never be polarized in air!

During this time, it is recommended to keep the protection cap filled with fresh conditioning solution on the sensor (see “Section 7 – Storage”), especially if the sensor is being polarized for longer than 6 hours (Exemption: the ISM version InPro 6850i reaches polarization after 1h and is then ready for operation).

A shorter polarization period is sufficient if the sensor has been disconnected for only a few minutes. The following table serves to establish the correct polarization time in relation to the depolarization time.

| Depolarization time ¹ t_{depol} [Min.] | Minimum required polarization time ² [Min.] | |
|---|---|--------------|
| | InPro 6900 i; InPro 6950 i | InPro 6850 i |
| > 30 | 360 | 60 |
| 15 to 30 | $6 \times t_{\text{depol}}$ | 45 |
| 5 to 15 | $4 \times t_{\text{depol}}$ | 30 |
| < 5 | $6 \times t_{\text{depol}}$ | 0 |

- 1 Depolarization time: Time span in which the polarization voltage is cut off from the sensor. This is the case:
 - during the time the cable is disconnected or no transmitter or sensor master is connected to the cable, or the transmitter has been disconnected from the current supply.
 - after changing the electrolyte or membrane body. In this case at least 6 hrs (InPro 6850i; 1 h) of polarization must follow.
- 2 Polarization time: Time span during which the sensor is under a polarization voltage.

Important! Setting of the polarization voltage on the transmitter for correct measurements:

- Standard applications InPro 6850i: **–675 mV**

- Measurement of permanently low oxygen concentrations (<500 ppb in liquids or <10,000 ppm [vol.] in gases) in the presence of volatile acidic components (e.g. carbon dioxide during measurements in breweries) e.g. InPro 6900 i/6950 i: **-500 mV**

Note: To ensure the supply of the correct polarization voltage the transmitter must be set accordingly.

5.2 Calibration

5.2.1 Purpose of calibration

Each oxygen sensor has its own individual slope and own individual zero point. Both values are subject to change, for example, through electrolyte consumption or after exchange of electrolyte or membrane body. To ensure high measurement accuracy of the sensor, a calibration must be carried out regularly, but at least after each change of electrolyte or membrane. Prior to calibration, the sensor has to be polarized for at least 6 hours (InPro 6850 i; 1 h).

Please remove the protection cap from the sensor, rinse the sensor with water and air dry it before starting calibration.

To check if your sensor needs a recalibration, you may dry it and take it in the air to check that the reading is close to 100%. If not, then the sensor needs a new calibration.

With ISM sensors all calibration data are stored in the sensor.

5.2.2 What you have to know for calibration

Calibration is generally carried out at the polarization voltage at which measurements are taken.

Zero point calibration is recommended:

- when measuring less than 5 ppb in liquids
- when measuring less than 125 ppm (by vol.) in gases

General remarks:

- For calibration in air, the sensor membrane must be dry, since adhering water drops can falsify the measured oxygen value.
- Make sure that the oxygen saturation index of the calibration medium is correct and remains constant during calibration.
- In the event of calibration in water or sample medium, the calibration medium must be in equilibrium with the air. Oxygen exchange between water and air is very slow. Therefore it takes quite a long time until water is saturated with atmospheric oxygen.

- For correct calibration, a minimum flow rate of the calibration medium is necessary.
- Make sure that all other parameters, such as temperature and pressure, are constant.

For continuous applications, we recommend **periodic recalibration in line with your requirements on accuracy, the type of process in operation and your own experience**. The frequency of the need for re-calibration depends very much on the specific application, and therefore appropriate intervals cannot be exactly defined here.

5.2.3 Single point calibration

By carrying out a single point calibration, the factual slope of the sensor can be established. The calibration medium can be water with known oxygen saturation index (e.g. air-saturated water) or air with known water-vapor saturation (e.g. water-vapor saturated air).

After the sensor signal has stabilized, the complete measuring system can then be calibrated to the 100 % value of the desired measurable variable, e.g. 100 % air, 20.95 % O₂, or 8.26 ppm at 25 °C (77 °F) and normal pressure (see instruction manual for the transmitter).

5.2.4 Dual point calibration

By carrying out a dual point calibration both slope and zero point of the sensor can be established.

Important! In case of a dual point calibration, always start by the zero point calibration before calibrating the slope.

As a rule, the zero point should be adjusted to zero manually, or it is automatically performed by the transmitter (see instruction manual for the transmitter).

Zero point calibration is recommended:

- when measuring less than 5 ppb in liquids
- when measuring less than 125 ppm (by vol.) in gases

Attention! Incorrect zero point calibration is a frequent source of measurement error. For correct calibration, we recommend the use of carbon dioxide gas with a level of purity of at least 99.9995 %.

After the sensor signal has stabilized (after **6...12 hours**), the sensor can be calibrated through the relevant transmitter to the zero % value of the desired measurable variable, e.g. 0 % air, 0.0 % O₂, or 0.0 ppm (see instruction manual for the transmitter).

6 Maintenance

6.1 Inspection of the sensor

6.1.1 Visual inspection

To check your sensor, we recommend the following procedure:

- The contacts of the connector must be dry. Moisture, corrosion and dirt in the connector can lead to false readings.
- Check the cable for buckling, brittle areas or ruptures.
- Before calibration always examine the membrane foil optically for signs of damage. The foil must be intact and clean. Dirty membranes should be wiped clean using a soft, moist tissue.

Note: An undulated membrane has no influence on the sensor performance, assuming the membrane is intact.

- The membrane body must be replaced if the sensor has too long a response time, the reading is unstable or subject to drift, and if the sensor cannot be calibrated or the membrane shows sign of mechanical damage.
- Check the cathode area for discoloration, contamination or cracks in the glass. If necessary rinse with demineralized water and clean with a clean soft brush or soft paper tissue.

Attention! Do not use any cleaning agents containing alcohol. This could damage the sensor or lead to a faulty current.


Attention! The glass body is fragile and sensitive to vibration.

6.1.2 Testing the METTLER TOLEDO O₂ sensor master

We recommend using the METTLER TOLEDO O₂ sensor master to check the quality of your sensor as follows:

- **Connect the sensor to the O₂ sensor master.**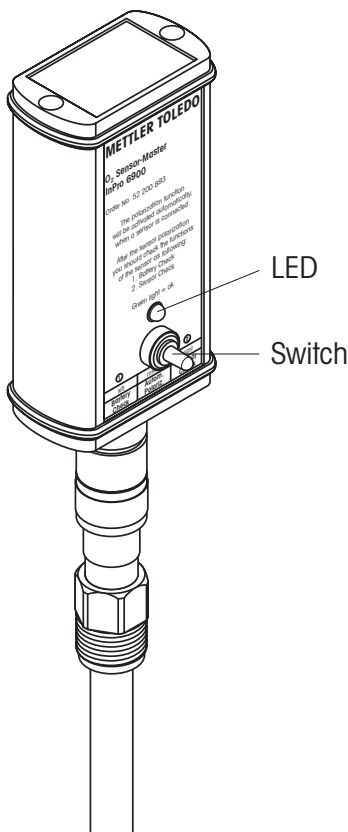

As soon as the sensor is connected to the O₂ sensor master, the polarization function is automatically activated. Please note: if the sensor was disconnected from the transmitter for longer than 5 minutes, the sensor must be polarized first (polarizing time see "Section 5.1") to get representative test results.

- **Battery Check:**
Push the switch to the left. If the battery is ok and the O₂ sensor master is operational the green LED lights up. Otherwise, please consult the instruction manual of the O₂ sensor master.
- **Sensor Check:**
For this test the O₂ sensor must be fully polarized and the membrane of the sensor must be dry and clean.

Expose the sensor (connected to O₂ sensor master) to the air. By pushing the switch to the right to the position "2 – Sensor check", the O₂ sensor master checks whether the electrode current for air measurement delivered by the sensor is within the admissible range, i.e. 2500 to 6000 nA for InPro 6950.

If the green LED lights up the current for air measurement is within the admissible range.

If the LED does not light up, you should check the battery of the O₂ sensor master (see instruction manual "Accessories"). If the battery is working, then there is probably a problem with the sensor. You should change the electrolyte and/or the membrane body of the sensor. If after a membrane change the LED still does not light up, this means that there is possibly something wrong with the interior body of the sensor. You should then change it (see "Section 6.2").

Important! The Sensor Check function only verifies the correctness of the electrode current for air measurement. In order to be absolutely sure of the functionality of the sensor, the residual signal in an oxygen free medium should also be controlled (see "Section 6.1.3").

6.1.3 Testing the sensor via a transmitter

A periodic zero current measurement is recommended for verification of proper sensor function.

Note: At the time you carry out the zero current measurement, the sensor must be polarized.

Zero current measurement can be done by using zeroing gel (order no. 30 300 435) or nitrogen (N₂) or carbon dioxide (CO₂) calibration gases with a purity of at least 99.995 %, or alternatively in a sample medium saturated with one of these gases.

After 2 minutes in an oxygen-free sample medium, the reading on the transmitter should drop to below 10 % of the reading in ambient air, and within 10 minutes the value should have dropped to below 1 %.

If the measured values are too high, this suggests a depleted electrolyte or a defective membrane. In the first instance replace the electrolyte, and in the second case exchange both the membrane body and the electrolyte accordingly.

If after such procedures the above mentioned values are still not reached, replace the interior body. If this does not solve the problem send the sensor to your local METTLER TOLEDO representative for inspection.

Many sample media contain volatile substances which, even at very low concentrations, have a clearly perceptible smell. Similarly to oxygen, these substances are able to invade the electrolyte through the gas-permeable membrane. Accordingly, they become noticeable when changing the electrolyte. In most cases, such substances have absolutely no influence on the measuring properties of the sensor. Slight discoloration of the electrolyte also has no effect on the measuring properties.

6.1.4 ISM design

The integrated ISM functionality allows extensive monitoring of the sensor. The following parameters are stored in the sensor:

- serial no.
- type of sensor
- order no.
- calibration data
- CIP/SIP counter
- slope
- zero point

6.2 Changing the electrolyte, the membrane body or the interior body

Note: The InPro 6900 i and 6950 i use a special electrolyte which contains an oxygen scavenger. This electrolyte ensures fast response time and together with the guard ring increases the signal stability of the sensor. The electrolyte must be exchanged on a regular basis or when the sensor has been exposed to the air for more than 24 hours without a watering cap filled with conditioning solution.

Note: The content of the electrolyte bottle has to be used within 3 months after opening.

If the membrane and/or the interior body exhibits signs of failure (long response time, increased current in an oxygen-free medium, mechanical damage, etc.) the membrane body and/or the interior body has to be replaced.

Warning! The O₂ electrolyte has an alkaline pH value. Contact of electrolyte with mucous membranes or eyes is to be avoided. Therefore protective gloves and safety glasses have to be worn for the following dismantling works. If such contact occurs, the affected area should be well rinsed with water. In the case of accident, or should ever any adverse signs appear, get immediate medical attention.

When changing the electrolyte, the membrane body or the interior body, please observe the following instructions (see also the following illustration):

Attention! Make sure that this maintenance step is carried out in a clean place.

1. Unscrew the cap sleeve from the sensor shaft and carefully pull it off the sensor.
2. Pull off the membrane body from the interior body. If it is tightly fitted, eject by pushing it with a flat finger tip. Before electrolyte is refilled, the membrane body must be removed from the cap sleeve!
3. Rinse the interior body with demineralized water and carefully dab it dry with a paper tissue.

Note: steps 4 to 7 may only be carried out when changing the interior body.

4. Unscrew the retainer nut of the interior body with an adjustable wrench or with a 9 mm wrench.
5. Remove the interior body by pulling it out of the sensor shaft. If necessary use pliers.


Warning! Do not twist the interior body. Otherwise the connection pins can be damaged.

6. Insert the new interior body in the sensor shaft. Turn the interior body in the shaft until the slit of the interior body is aligned with the pin placed in the shaft.
7. Press the body in the shaft and screw the new retainer nut in place.
8. Examine the O-rings visually for mechanical defects, and replace if necessary.
9. Half-fill the new membrane body with O₂ electrolyte.

Note: The electrolyte bottle is equipped with a special pouring system. To ensure proper functioning, hold the bottle vertically, upside-down.

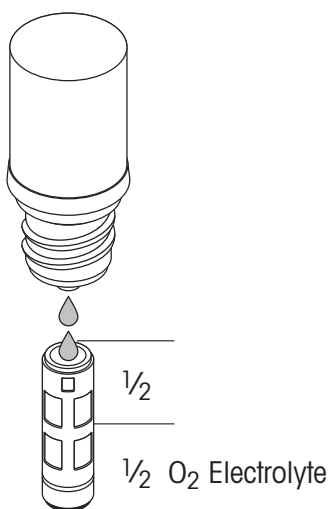Note: Make sure that all air bubbles are removed from the membrane body. Air bubbles can be removed by carefully tapping on the membrane body.


10. Slip the membrane body over the interior body while holding the sensor in a vertical position. Excess electrolyte will be displaced and has to be removed with a paper tissue.

Important! No electrolyte, sample media or contamination may be present between the membrane body and the cap sleeve. Please check carefully!

11. Carefully slip the cap sleeve over the fitted membrane body, holding the sensor in a vertical position and screw it tight. The cap sleeve must be clean and dry.
12. After each exchange of electrolyte or membrane body, the sensor has to be repolarized and recalibrated.

Replacement of the interior body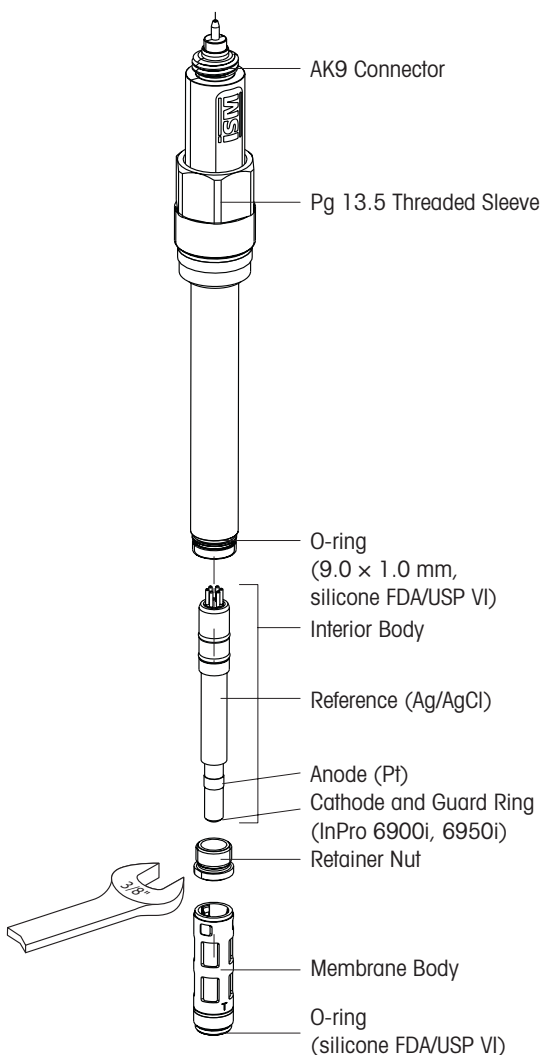

7 Storage

For storage periods longer than 24 hours, we recommend you use a protection cap filled with the same cleaning and conditioning solution as that used for our portable DO measurement system (order no. 52 200 255). This solution contains an oxygen scavenger which ensures that the electrolyte of the sensor does not come into contact with oxygen when not in use.

When the sensor is stored without polarization for more than one week, the electrolyte has to be removed.

Prepare the cleaning and conditioning solution as follows: Dissolve one tablet in 40 ml of deionized water and wait 5 minutes for the tablet to be completely dissolved. Fill the protection cap with this solution and place it over the tip of the sensor. This solution has some cleaning properties which will keep the membrane free of microorganisms. In case you do not have a cleaning and conditioning set, you may also use zeroing gel or deaerated water in the protection cap. Before mounting the sensor in the process, always remove the protection cap and rinse the tip of the sensor with water.

Attention! If the storage period of the sensor without current supply (transmitter, sensor master) exceeds 1 week, the sensor should be stored dry, i.e. without any electrolyte in the membrane body. A sensor being stored dry (without electrolyte in the membrane body) may on no account be connected to the O₂ sensor master or any other polarization module.

8 Product specification

8.1 Certificates

Each sensor is delivered with a set of **3.1 certificates** (complying with EN 10204).

All wetted metal parts (sensor shaft, cap sleeve and membrane body) are identified with an engraved symbol corresponding to the heat number on the paper certificate delivered with the sensor.

Each wetted metal part (sensor shaft, cap sleeve and membrane body) is polished in order to get a surface roughness lower than 0.4 µm (16 µin). This represents a roughness grade number of N5 (according to ISO 1320:1992).

8.2 Specifications

| InPro 6850 i/6900 i/6950 i | |
|---|---|
| Measurement principle | Polarographic Clark electrode |
| Working conditions | |
| Pressure resistance (measurement) | 6850i: 0.2 ... 6 bar 6900i: 0.2 ... 6 bar (9 bar with T-6900R) 6950i: 0.2 ... 6 bar |
| Mechanical pressure resistance | Max. 12 bar |
| Temperature range (measurement) | 0 ... 80 °C (32 ... 176 °F) |
| Temperature range (mechanical) | 6850i: - 5 ... 140 °C (23 ... 284 °F) 6900i: - 5 ... 121 °C (23 ... 249.8 °F) [sterilizable and autoclavable] 6950i: - 5 ... 121 °C (23 ... 249.8 °F) [sterilizable] |
| Construction | |
| Temperature compensation | Automatic |
| Cable connection | AK9 |
| O-ring material | Silicone FDA and USP Class VI approved |
| Membrane material | PTFE/Silicone/PTFE (reinforced with steel mesh) |
| Material sensor body (wetted parts) | 316L stainless steel |
| Surface roughness of wetted metal parts | N5 (R _a = 0.4 µm [16 µin]) |
| Quick disconnect interior body | Standard |
| Cathode | Pt |
| Anode | 6850i: Pt 6900i: Ag 6950i: Pt |
| Guard ring | 6850i: - 6900i: Pt 6950i: Pt |
| Reference | Ag |
| Dimensions | |
| Sensor diameter | 6850i: 12/25 mm 6900i: 12 mm 6950i: 12 mm |
| Immersion length (a) for 12 mm sensor | 6850i/6900i: 70, 120, 220, 320, 420 mm (2.8, 4.7, 8.66, 12.6, 16.5") 6950i: 70, 120, 220, 320 mm (2.8, 4.7, 8.66, 12.6") |
| Immersion length (a) for 25 mm sensor | 6850i: 80, 160, 260, 360 mm (3.2, 6.3, 10.2, 14.2") 6900i: - 6950i: - |
| Performance | |
| Detection limit | 6850i: 6 ppb to saturation 6900i: 1 ppb to saturation 6950i: 0.1 ppb to saturation |
| Accuracy | 6850i: ≤ ± (1 % + 6 ppb) 6900i: ≤ ± (1 % + 1 ppb) 6950i: ≤ ± (1 % + 0.1 ppb) |
| Response time at 25 °C/77 °F | 98 % of final value in < 90 s |

| | |
|---|--|
| Sensor signal in ambient air at 25 °C/77 °F | 6850 i: 50 ... 110 nA 6900 i: 250 ... 500 nA 6950 i: 2500 ... 5000 nA |
| Residual signal in oxygen-free medium | 6850 i: <0.1 % of the signal 6900 i: <0.3 % of the signal 6950 i: <0.025 % of the signal |

Certification

| | |
|----------------------------------|-----------------------------------|
| EHEDG, 3A | Yes |
| 3.1 (EN 10204) | Yes |
| ATEX, IECEx and UKEX Certificate | 6850 i/6900 i: Yes 6950 i: Yes |
| FM Approval | 6850 i/6900 i: Yes 6950 i: Yes |
| FDA/USP VI | Yes |
| Quality Control | Yes |

Compatibility

| | |
|----------------------------------|-------------------|
| with METTLER TOLEDO transmitters | see "Section 9.4" |
| with METTLER TOLEDO housings | see "Section 9.5" |

9 Ordering information

For more detailed information refer to the technical data sheet or ask your local distributor.

9.1 Sensors with ISM functionality

Ordering Information

| – straight version | 6850i | 6900i | 6950i |
|--------------------|------------|------------|------------|
| 70 mm | 52 206 118 | 52 206 316 | 52 206 127 |
| 120 mm | 52 206 119 | 52 206 317 | 52 206 128 |
| 220 mm | 52 206 120 | 52 206 318 | 52 206 129 |
| 320 mm | 52 206 121 | 52 206 319 | 52 206 130 |
| 420 mm | 52 206 122 | – | – |
| – 25 mm version | | | |
| 80 mm | 52 206 123 | – | – |
| 160 mm | 52 206 124 | – | – |
| 260 mm | 52 206 125 | – | – |
| 360 mm | 52 206 126 | – | – |

9.2 Accessories

| Accessories | Order No. |
|--|------------|
| O ₂ sensor master digital ISM | 52 206 329 |
| Cables | |
| AK9 coax cable with K8S connector 1 m | 59 902 167 |
| AK9 coax cable with K8S connector 3 m | 59 902 193 |
| AK9 coax cable with K8S connector 5 m | 59 902 213 |
| AK9 coax cable with K8S connector 10 m | 59 902 230 |
| AK9 coax cable with K8S connector 20 m | 52 300 204 |

9.3 Spare parts

| Spare parts | Order No. |
|--|------------|
| – InPro 6850i | |
| Membrane body (single), T-96 | 52 200 071 |
| Membrane kit T-96 (4 membranes, 1 O-ring set silicone, 25 ml of electrolyte, wetted parts SS 316L) | 52 200 024 |
| Membrane bodies (16 pcs), T-96 | 52 206 114 |
| O ₂ Electrolyte pack (3 × 25 ml) | 30 298 424 |
| Interior body (replacement), InPro 6850i | 52 206 347 |
| – InPro 6900i | |
| Reinforced membrane body (single), InPro 6900 (T-6900R) | 52 201 108 |
| Reinforced membrane kit InPro 6900 (T-6900 R) | 52 201 109 |
| InPro 6900 electrolyte pack (3 × 5 ml) | 30 298 425 |
| Interior body (replacement), InPro 6900 | 52 200 943 |
| – InPro 6950i | |
| Membrane body (single) | 52 206 105 |
| Membrane kit (4 membranes, 1 O-ring set, 2 × 5 ml of electrolyte) | 52 206 106 |
| InPro 6950 electrolyte pack (3 × 5 ml) | 30 298 426 |
| Interior body (replacement), InPro 6950 | 52 206 112 |

9.4 Recommended transmitters

| Transmitters M700 | Order No. |
|--|------------------|
| Transmitter base, coated, M700C | 52 121 171 |
| Transmitter base, coated, Ex, VPW, M700XC/VPW | 52 121 172 |
| Transmitter base, coated, Ex, 24V, M700XC/24V | 52 121 173 |
| Transmitter base, stainless steel, M700S | 52 121 174 |
| Transmitter base, stainless steel, Ex, VPW, M700XS/VPW | 52 121 175 |
| Transmitter base, stainless steel, Ex, 24V, M700XS/24V | 52 121 176 |

| Transmitter modules M700 | Order No. |
|--|------------------|
| ISM O ₂ module 4700i | 52 121 263 |
| ISM O ₂ module, Ex, 4700iX | 52 121 263 |
| ISM O ₂ module 4700i ppb | 52 121 265 |
| ISM O ₂ module, Ex, 4700iX ppb | 52 121 266 |
| ISM O ₂ module 4700i traces | 52 121 295 |
| ISM O ₂ module, Ex, 4700iX traces | 52 121 294 |

| Transmitters M300D | Order No. |
|--|------------------|
| M300i, ¼DIN, single-channel, multi-parameter | 52 121 354 |
| M300i, ½DIN, single-channel, multi-parameter | 52 121 355 |
| M300i, ¼DIN, dual-channel, multi-parameter | 52 121 356 |
| M300i, ½DIN, dual-channel, multi-parameter | 52 121 357 |

| Transmitters M400 | Order No. |
|--------------------------|------------------|
| M400, type 1 | 52 121 348 |
| M400, type 2 | 52 121 349 |
| M400, type 3 | 52 121 350 |

For more information consult the "Parameter Fit Guide" below.

Parameter Fit Guide for Transmitter M400

| | M400 | M400 | M400 |
|-----------------------------|---------------|---------------|---------------|
| Conventional sensors | Type 1 | Type 2 | Type 3 |
| pH/ORP | • | • | • |
| Cond 2-electrode | • | • | • |
| Cond 4-electrode | • | • | • |
| DO ppm | – | • | • |
| DO ppb | – | – | • |
| ISM sensors | | | |
| pH/ORP | • | • | • |
| Cond 2-electrode | • | • | • |
| Cond 4-electrode | • | • | • |
| DO ppm | – | • | • |
| DO ppm optical | – | – | • |
| DO ppb | – | – | • |
| DO trace | – | – | • |

9.5 Recommended housings

Housing (12 mm Ø)

Static housing

InFit 761 CIP

Retractable housings

InTrac 777 e

InTrac 797 e

Immersion housing

InDip™ 550

Note: The housings are available in different versions. Please contact your local METTLER TOLEDO sales organization to get the correct ordering information.

10 Theory of the polarographic sensor

10.1 Introduction

Two types of electrodes are employed in analytical sensors **potentiometric** and **amperometric** electrodes.


- **Potentiometric electrodes** develop a voltage generated by the activity of a particular ion. Examples of such electrodes are glass electrodes (like pH electrodes) and most ion-selective electrodes. Their individual potentials cannot be determined. The measurable quantity is the difference of potential between the measuring electrode and an inert reference electrode. The potential of the reference electrode must be constant.

All potentiometric electrodes are subject to Nernst's law and for this reason electrodes and measuring instruments are in most cases interchangeable. An important requirement of potentiometric measurements is the virtually currentless determination of the electrode voltage. During measurement no chemical reaction occurs and the solution remains in equilibrium.

- In the case of **amperometric electrodes**, such as the **oxygen electrode**, activity measurement is based on a current measurement.

The conventional oxygen electrode consists of a cathode and an anode conductively connected by an electrolyte. A suitable polarization voltage between the anode and the cathode selectively reduces the oxygen at the cathode.

Cathode reaction: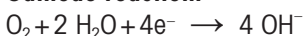

These chemical reactions result in an electric current which is proportional to the oxygen partial pressure ($p\text{O}_2$). The oxygen electrode constantly reduces oxygen. Thereby, the concentration of dissolved oxygen is reduced. By diffusion this oxygen is replaced. The viscosity and flow rate of the solution are therefore important parameters. The electrode current of an oxygen electrode is determined not only by the oxygen partial pressure but by many other electrode parameters. The electrode currents of different electrode types may differ by several powers of ten. For this reason oxygen electrodes and amplifiers cannot be freely interchanged.

10.2 Principle of the design of an oxygen electrode

There are two main types of oxygen electrode:

- Electrodes **without** membrane
- Electrodes **with** gas-permeable membrane (Clark Principle)

The membrane electrode according to Clark is today most widely used. As compared to the electrode without membrane it possesses the following advantages:

- Oxygen measurement in gases and solutions
- No mutual contamination of electrode and solution
- No or little dependency on flow



In the case of the Clark electrode, geometrical configuration is very important. In particular, the thickness of the electrolyte film between the cathode and the membrane must be within narrow tolerances so as to ensure good linearity and a low zero current (current in a nitrogen atmosphere).

METTLER TOLEDO oxygen sensors are available in different designs:

Type A, 2 electrode system, InPro 6800

InPro 6800 for medium and high oxygen concentrations. Cathode and anode/reference. Anode and reference are united in a silver/silver chloride electrode.

The following equilibration reaction takes place at the anode resp. reference: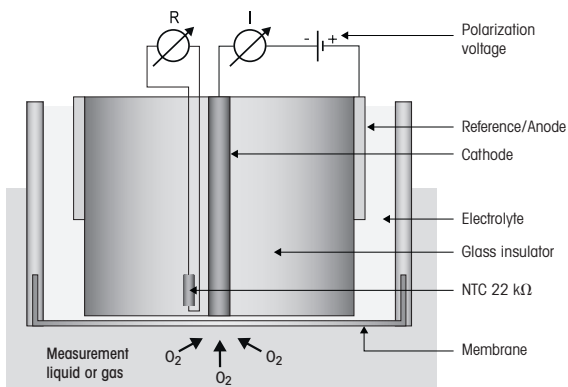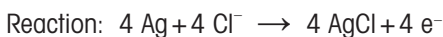

Type B, 3 electrode system, InPro 6850 i

InPro 6850i for medium and high oxygen concentrations. The reference corresponds the conventional silver/silver chloride anode. The anode is a platine electrode and separated from the reference.


The following reaction takes place at the anode:

The reference is a silver/silver chloride electrode. After polarization there is a stable equilibrium between electrode surface and electrolyte. No net-reaction takes place.


Reference reaction in equilibrium:

Reaction: No current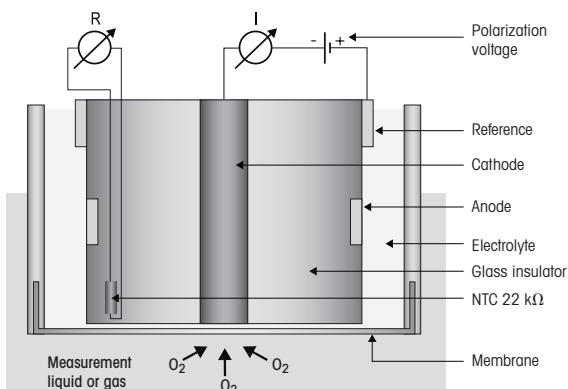



Type C, 3 electrode system, InPro 6900 (i)

InPro 6900 and InPro 6900 i. For oxygen measurements in the lower ppb range. Here, the anode and reference are united in a silver/silver chloride electrode (as with type A). The sensors are equipped with an additional guard ring around the cathode. Like the cathode with the anode, this forms a closed electric circuit, which prevents oxygen diffusing from the side to the cathode and falsifying the measurement result.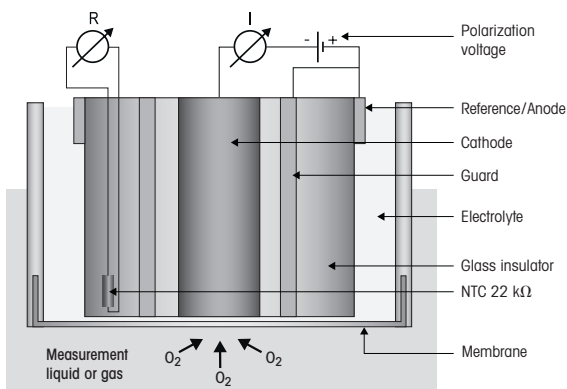


In all sensors the function of the cathode is identical.

Between cathode and reference a constant voltage of -500 resp. -675 mV (polarization voltage) is applied. Thereby a reduction of O_2 into OH^- at the cathode takes place.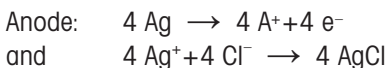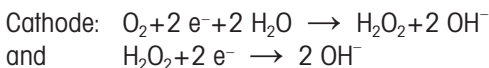



The total equation thus is:

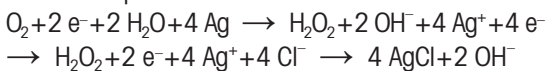
Type D, 4 electrode system, InPro 6950 (i)

The InPro 6950(i) is intended for measuring permanent low oxygen levels down to 0.1 ppb. In terms of measurement principle it is a combination of the InPro 6850i and InPro 6900(i). It consists of 4 electrodes. The anode (platinum) and reference (silver/silver chloride) are separated in two electrodes. A guard ring is placed around the cathode. The cathode has the highest diameter of all amperometric oxygen sensors from METTLER TOLEDO.


The electrodes show the following reactions: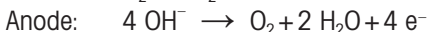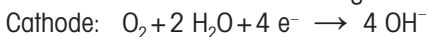

Reference: No current

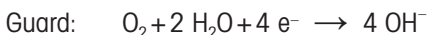Due to the larger surface of the cathode a higher current appears during measurement. This allows for measurement of oxygen levels below 1 ppb.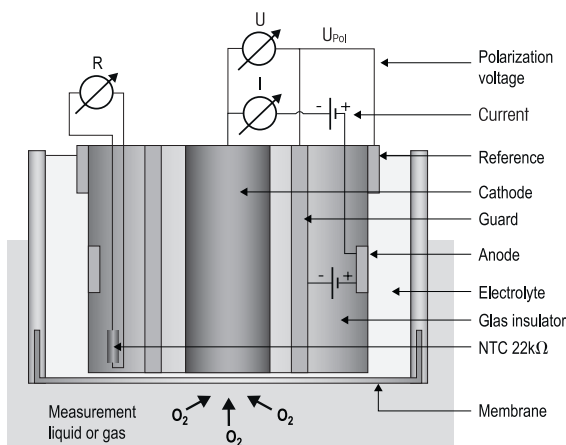

As a consequence of the reaction at the cathode and the anode, a current flows between cathode and anode (4 electrons per oxygen molecule), which is proportional to the oxygen partial pressure at the cathode.

The amperage of the current flow depends on the surface area of the cathode. Typical values are:

| | Air current | Zero current in % of the air current |
|----------|--------------------|---|
| 6800 | 50–110 nA | <0.1 |
| 6850 (i) | 50–110 nA | <0.1 |
| 6900 (i) | 250–500 nA | <0.03 |
| 6950 i | 2500–5000 nA | <0.025 |

This current is measured and converted and indicated in the transmitter into an oxygen value. In ISM sensors this computation takes place in the sensor and the transmitters displays this value.

10.3 Parameters determining current

The quantity of oxygen diffused in and the magnitude of the electrode current are influenced by the following parameters:

- Oxygen partial pressure of the solution
- Membrane material and thickness
- Size of cathode
- Polarization voltage
- Temperature
- Flow conditions in the solution


Fick's law gives the mathematical relationship between these parameters:

$$I = k \times D \times a \times A \times \frac{pO_2}{x}$$

- I = Electrode current
- k = Constant
- D = Coefficient of diffusion of O₂ in the membrane
- a = Solubility of O₂ in the membrane material
- A = Cathode surface
- pO₂ = Oxygen partial pressure in the solution
- x = Thickness of gas-permeable membrane

10.4 Polarization voltage

The voltage between the anode and the cathode is so selected that the oxygen is fully (> A, see polarogram) reduced while other gases are unaffected (< D). The ideal voltage for the Pt/Ag/AgCl or Pt/Pt/Ag/AgCl system is between –500 and –750 mV.

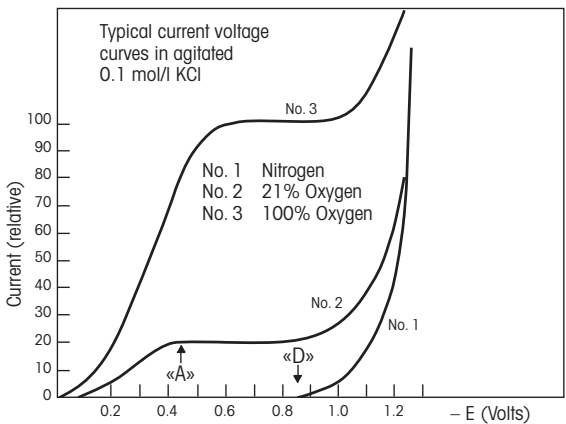The polarization voltage should remain as constant as possible. Besides a constant source of voltage the following prerequisites must be met: The electrical resistance of electrolyte film must not exceed a particular value so as to avoid a voltage drop.

The anode must have a large surface so as to prevent polarization of the anode by the electrode current.

10.5 Temperature

The temperature dependence of the current passing through an oxygen electrode when referred to a constant O₂ partial pressure is determined mainly by the properties of the gas-permeable membrane.

10.6 Dependence on flow

With most oxygen electrodes the electrode current is smaller in stagnant solutions than in agitated ones. In consequence of the oxygen consumption of the electrode, a reduction of oxygen takes place outside the membrane in close proximity to the cathode. The missing oxygen is replaced by diffusion. If the electrode current is strong, the solution cannot fully restore the reduced oxygen by diffusion. This results in an electrode current weaker than would correspond to conditions in the solution. In agitated solutions the oxygen is conveyed to the surface of the membrane not only by diffusion but additionally by the flow (convection). In that case no oxygen impoverishment occurs at the membrane surface.

A high degree of flow dependence occurs mainly with large cathodes, and thin and highly permeable membranes, i.e. where electrode currents are large.

The problem of flow dependence is often solved by prescribing a minimum flow rate.

In METTLER TOLEDO InPro 6950 sensors, the thin PTFE membrane determining the electrode current (i.e. the actual measuring signal) is separated from the sample solution by a relatively thick silicone membrane. The latter is highly permeable to oxygen molecules and thus acts as an oxygen reservoir. The diffusion of oxygen out of the sample solution into the silicone membrane is spread over a wide area. Since this results in less oxygen being extracted from the sample solution per unit area, the double PTFE/silicone membrane forms an effective buffer against disturbances due to hydrodynamic flow.

This membrane coupled with the guard ring and the special electrolyte ensures excellent signal stability even when the hydrodynamic flow stops (as on a beer filling line for example).

10.7 Oxygen partial pressure – oxygen concentration

The electrode current is dependent on the partial pressure of oxygen and the oxygen permeability of the membrane. Conversion of partial pressure into concentration of oxygen depends on the measurement medium (measurement in liquids or gases).

Measurement in liquids

When measuring in liquids, the concentration of oxygen also depends on the solubility of the oxygen

in the measurement medium. However, since this is not measured by the sensor current, the concentration of oxygen must be calculated in the transmitter. To do this, Henry's law is applied which states that the concentration of oxygen is proportional to the partial pressure of oxygen (pO_2).

$$CI = pO_2 \times \alpha$$

α = Solubility factor

If " α " is constant, the oxygen concentration can be determined by means of the electrode. This applies at constant temperature and with dilute aqueous solutions such as drinking water.

The solubility factor is strongly influenced not only by the temperature but also by the composition of the solution:

| Medium, saturated with air | Solubility at 20 °C (68 °F) and 760 mm Hg |
|---------------------------------------|--|
| Water | 9.2 mg O ₂ /l |
| 4 mol/l KCl | 2 mg O ₂ /l |
| 50 % Methanol-water | 21.9 mg O ₂ /l |

Although the solubilities vary widely, the oxygen electrode gives the same reading in all three solutions.

Thus, determination of the oxygen concentration is only possible with constant and known solubility factors " α ".

Solubility may be determined by a Winkler titration or the method developed by Käppeli and Fiechter.

Measurement in gases

The concentration of oxygen when measuring in gases is always given as a proportion by volume of the gas composition. Common units are % (by vol.) ppm (by vol.).

They can simply be converted from one unit of measurement to the other.

Example:

The percent by volume of the composition of air is generally known. For example, air contains 20.95 % oxygen. This corresponds to 209,500 ppm (by vol.). (Conversion: ppm value = 10,000 × value in %)

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US-Normen

| Titel | Nummer | Ausgabe- datum |
|--|--------------------------|---------------------------|
| Approval Standard for Electrical Equipment for Use in Hazardous (Classified) Locations – General Requirements | FM Class 3600 | 2022 |
| Approval Standard for Intrinsically Safe Apparatus and Associated Apparatus for Use in Class I, II & III, Division 1, Hazardous (Classified) Locations | FM Class 3610 | 2021 |
| Approval Standard for Electrical Equipment for Measurement, Control and Laboratory Use | FM Class 3810 | 2021 |
| Explosive Atmospheres – Part 0: Equipment – General Requirements | ANSI/UL 60079-0:2020 | 2020 |
| Explosive Atmospheres – Part 11: Equipment Protection by Intrinsic Safety “i” | ANSI/UL 60079-11:2018 | 2018 |
| Safety Requirements for Electrical Equipment for Measurement, Control, and Laboratory Use – Part 1: General Requirements | ANSI/UL 61010-1:2018 | 2018 |

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| Titel | Nummer | Ausgabe- datum |
|--|--|---------------------------|
| Explosive Atmospheres – Part 0: Equipment – General Requirements | CAN/CSA-C22.2 No. 60079-0:2019 | 2019 |
| Explosive Atmospheres – Part 11: Equipment Protection by Intrinsic Safety “i” | CAN/CSA-C22.2 No. 60079-11- 14:R2018 | R2018 |
| Safety Requirements for Electrical Equipment for Measurement, Control, and Laboratory Use – Part 1: General Requirements | CAN/CSA-C22.2 No. 61010-1- 12:R2017 | R2017 |

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| Titre | Numéro | Date de publication |
|--|-----------------------|----------------------------|
| Approval Standard for Electrical Equipment for Use in Hazardous (Classified) Locations – General Requirements | FM Class 3600 | 2022 |
| Approval Standard for Intrinsically Safe Apparatus and Associated Apparatus for Use in Class I, II & III, Division 1, Hazardous (Classified) Locations | FM Class 3610 | 2021 |
| Approval Standard for Electrical Equipment for Measurement, Control and Laboratory Use | FM Class 3810 | 2021 |
| Explosive Atmospheres – Part 0: Equipment – General Requirements | ANSI/UL 60079-0:2020 | 2020 |
| Explosive Atmospheres – Part 11: Equipment Protection by Intrinsic Safety “i” | ANSI/UL 60079-11:2018 | 2018 |
| Safety Requirements for Electrical Equipment for Measurement, Control, and Laboratory Use – Part 1: General Requirements | ANSI/UL 61010-1:2018 | 2018 |

Normes canadiennes

| Titre | Numéro | Date de publication |
|--|-------------------------------------|----------------------------|
| Explosive Atmospheres – Part 0: Equipment – General Requirements | CAN/CSA-C22.2 No. 60079-0:2019 | 2019 |
| Explosive Atmospheres – Part 11: Equipment Protection by Intrinsic Safety “i” | CAN/CSA-C22.2 No. 60079-11-14:R2018 | R2018 |
| Safety Requirements for Electrical Equipment for Measurement, Control, and Laboratory Use – Part 1: General Requirements | CAN/CSA-C22.2 No. 61010-1-12:R2017 | R2017 |

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All rights reserved.